



## Shoreline Complete™ V4 Protocol 2

Lyse, Purify, & Amplify Sets C1, C2, C3, C4, C5, C6 (16 Tubes)  
Saliva, Swabs, and Cell Pellets

Rapidly purify microbial DNA from low input samples (e.g., saliva, skin swabs, cell pellets) and prepare for sequencing using single step PCR amplification of the of the 254-base pair V4 region of the 16S rRNA gene for microbial community profiling. This sample kit is one of six that contain unique barcoded primer sets for processing 16 samples, each primer has the structure: **<TruSeq Universal Adapter: Part A><8 base barcode><TruSeq Universal Adapter Part B><16S specific primer>**. Every sample will have a unique barcode at the completion of the PCR reaction for sequencing all the samples together.

### Products

1. Shoreline Complete™ V4 Lyse, Purify, & Amplify Set C1...C6 (16 tubes) Cat. No. SCV4-C1...C6

### Additional Items Needed for 16 Barcode Kits: (Not included; see Suggested Supplier list below)

1. Additional Items Required for Lyse and Purify
  - a. 8-strip 0.2 mL PCR tubes w/individual domed caps for sample lysis and purification
  - b. Sterile Laboratory Grade Water
  - c. Fresh 70% ethanol
  - d. TE Buffer - 10mM Tris, 1mM EDTA, pH 8.0
  - e. 0.4M KOH Solution
  - f. Magnetic Rack ('bar' or 'post' suggested for higher DNA yield)
2. Additional Items Required for Amplify:
  - a. Qiagen MinElute® PCR Purification Kit
  - b. Sterile Laboratory Grade Water
  - c. Sterile 100% Ethanol
3. Suggested items for rapid processing
  - a. Repeater Pipette with multi-volume tips
  - b. Single Channel Manual Pipettes
  - c. Multichannel Pipette 0.5-10 µL working volume
  - d. Centrifuge with 96-well plate adaptor
  - e. FLOQSwabs Sterile Collection Swabs 30mm (16)
4. Suggested items for sample QC
  - a. 0.8% agarose gel with ethidium bromide or other dsDNA stain
  - b. Gel loading dye to mix with samples (e.g., NEB B7024)
  - c. 2kb DNA ladder

### Kit Contents for Shoreline Complete™ Lyse and Purify Protocol

1. Two 8-tube strips of dried Lysis-1
2. Purification Buffer (two 1.5 mL tubes with **green cap**)

### Kit Contents for Shoreline Complete™ Amplify Protocol

1. 96-well PCR plate with 16 barcoded primers in wells A1-H2 (Columns 1 and 2)
2. 2X PCR Premix (**blue cap**)
3. PCR sealing film (1 sheet)



**Important Considerations for Lyse and Purify:**

1. Samples should be thawed and stored on ice.
2. Shoreline Biome **lysis is pH dependent**. Samples stored in buffer with pH < 6.5 should use **at most 10** µL input volume, otherwise the risk of suboptimal lysis affecting downstream results is high.
3. Samples may contain pathogens. Handle samples using Universal Precautions. Please wear personal protective equipment and observe standard laboratory practices for handling, cleaning, and disposing of materials that come in contact with sample.
4. Shoreline Biome recommends including positive and negative controls in your project to monitor lysis and amplification efficiency as well as environmental contamination. Consider using mock microbiome cells (input=10<sup>9</sup>) and mock microbiome DNA (input 10-100ng) as positive controls and water as negative control.
5. Shoreline Biome lysis results in ssDNA; for quantitation, please choose an appropriate kit (e.g., Qubit ssDNA assay kit #Q12012, Promega Quantifluor ssDNA system, #E3190)

**Shoreline Complete™ Lyse and Purify Protocol**

1. **Make 0.4M KOH.** Measure 0.135g KOH into 15 mL tube, then add 6 mL H<sub>2</sub>O. 0.4M KOH expires after 7 days.
2. **Make 70% Ethanol.** Add 17.5 mL 100% ethanol to 7.5 mL sterile laboratory grade water
3. Remove **Purification Buffer (Green Cap)** from box and bring to room temperature.
4. **Sample Type:**
  - 4.1. **For Saliva:**
    - 4.1.1. Centrifuge saliva samples at 5500 RCF for 10 minutes to pellet cells.
    - 4.1.2. Remove and discard supernatant.
    - 4.1.3. Re-suspend in 50 µL of Molecular Biology Grade water, TE, or PBS.
    - 4.1.4. **Transfer 50 µL** of cell suspension (between 10<sup>6</sup> and 10<sup>8</sup> cells) to Lysis-1 tube. Do not mix.
  - 4.2. **For Skin swabs:**
    - 4.2.1. Moisten swab with buffer containing Tris-EDTA and 0.5% Tween 20.
    - 4.2.2. Collect sample from skin. Place swab in 1.7 mL microcentrifuge tube. Break the swab handle off at designated location.
    - 4.2.3. Centrifuge tube with swab for 5 minutes at 10000 RCF.
    - 4.2.4. Remove swab and transfer supernatant to a clean 1.7 mL tube.
    - 4.2.5. Centrifuge tube with swab second time for 5 minutes at 10000 RCF.
    - 4.2.6. Remove swab and combine supernatant with the first supernatant. Discard spent tube and swab.
    - 4.2.7. **Transfer 50 µL** of the supernatant to Lysis-1 tube. Do not mix.
  - 4.3. **For Pelleted cells:**
    - 4.3.1. Reconstitute the cell pellet by adding 50 µL Molecular Biology Grade water, TE, or PBS and pipetting up and down to re-suspend cells.
    - 4.3.2. **Transfer 50 µL of cell suspension** to Lysis-1 tube. Do not mix.
5. **Add 50 µL 0.4M KOH** to each tube of Lysis-1/sample mix. Do not mix. (A precipitate will form.)
6. Cap samples tightly; set PCR machine for proper tube height, close and lock the lid, and **heat samples to 95°C for 5 minutes**. Precipitate will dissolve. After incubation, **allow sample to cool to RT before unlocking PCR lid** – failure to do so may result in lids popping off tubes.
7. **While samples are cooling:**
  - 7.1. Label **clean 200 µL PCR tubes** for each sample

## Shoreline Complete™ V4 Protocol 2

Lyse, Purify, & Amplify Sets C1, C2, C3, C4, C5, C6 (16 tubes) Saliva, Swabs, and Cell Pellets

- 7.2. Vortex Purification Buffer tubes (**green cap**) to re-suspend any brown magnetic beads that may have settled
8. Add **100 µL of Purification Buffer (Green Cap)** to each sample. If precipitate formed, it should re-dissolve. Pipette up and down to mix.
9. Cap samples tightly; set PCR machine for proper tube height, close and lock the lid, and **heat samples to 50°C for 3 minutes** to allow DNA to bind to beads.
10. **After 50°C incubation, place samples on magnetic rack** to pellet beads with bound DNA to side of the tube (~60 seconds).
11. Once all beads are pelleted, **remove all supernatant SLOWLY and discard**. Avoid aspirating beads.
12. **Remove tubes from magnet, add 100 µL 70% ethanol** (at room temperature) to each tube, and re-suspend beads by pipetting gently up and down.
13. **Return tubes to magnet and allow beads to pellet** (~30 seconds)
14. **Pipette off residual ethanol and let air dry for 2 - 3 minutes**. Over-drying will make eluting DNA more difficult, so watch to ensure ethanol is mostly gone but bead pellet does not begin to shrink and crack.
15. **Remove tubes from magnet**.
16. Add **20 µL of 1X TE buffer** to tubes and briefly pipette up and down to re-suspend beads. Solution should turn brown in color to reflect beads mixed homogeneously throughout.
17. **Let solution incubate for 2 minutes at room temperature**.
18. **Place tubes on magnetic rack** for 30 - 60 seconds to **pellet magnetic beads. DNA is now in solution**.
19. Transfer supernatant containing eluted DNA into a **clean**, labeled microcentrifuge tubes on ice, avoiding pelleted beads.
20. **Remove tubes with pelleted beads from rack**.
21. **Add another 20 µL 1X TE buffer** onto bead pellets, gently pipette up and down to suspend beads.
22. **Incubate tubes for 2 minutes at room temperature**.
23. Return tubes to magnetic rack. Allow beads to pellet to sides for 30 - 60 seconds.
24. **Remove supernatant containing DNA** and combine with first eluate in the labeled tubes, avoiding pelleted beads. This is your **Shoreline Biome Lysis Mix**.

**NOTES:** DNA is **denatured**, due to KOH treatment during lysis. Although Lysis Mix may retain contaminants, the Shoreline Complete™ Amplify protocol is designed to work in the presence of these contaminants. Store Lysis Mix at +4°C for up to 10 days, or frozen at -20°C for 3 months. Avoid repeated freeze/thaws.

### Shoreline Complete™ Amplify Protocol

**NOTE: Only Columns 1 and 2 will be used.**

1. Remove the plate seal from the Amplicon Primer plate.
2. **Add 10 µL 2X PCR Premix (blue cap)** to Columns 1 and 2 of the 96-well plate.
3. **Add 10 µL Shoreline Biome Lysis Mix** to each of Column 1 and 2 wells.
4. Using the new PCR sealing film provided with the kit, cover the plate and seal tightly with no gaps.
5. Primers will dissolve off the bottom of each well. Observe that blue color is uniformly distributed throughout 20 µL reaction, if not, tap plate gently until uniform blue color is achieved in each well.
6. Spin plate briefly if reaction mix is not completely at the bottom of the wells.
7. Transfer the plate to PCR machine; close and lock the lid. Run PCR protocol:

## Shoreline Complete™ V4 Protocol 2

Lyse, Purify, & Amplify Sets C1, C2, C3, C4, C5, C6 (16 tubes) Saliva, Swabs, and Cell Pellets

95°C, 3:00	Denaturation
Repeat 34X	95°C, 0:20
	58°C for 0:15 ( <b>Ramp speed 4°C/sec</b> )
	72°C, 0:15
72°C, 2:00	Final extension

8. Check reactions on gel: Run 1.5 µL of sample with 5 µL diluted gel loading dye on 0.8% agarose gel in TBE, 150V for approximately 45 minutes, with DNA ladder. Band should be present at ~ 450bp.

### **Qiagen MinElute® PCR Purification Protocol**

1. Follow MinElute instructions for adding Ethanol (96% - 100%) to PE buffer, and label tubes.
2. Pool 5 µL of each amplified sample from the 96-well plate into a clean 1.5 mL microcentrifuge tube. (**Save 2 µL of pooled unpurified DNA for gel comparison with purified DNA in the last step below**)
3. Combine **78 µL** of pooled sample with 400 µL of Qiagen PB buffer.
4. Transfer **all** the resulting solution into the MinElute column with provided 2 mL collection tube.
5. Centrifuge the MinElute column at 17,900 RCF for 1 minute or until solution has passed through the column. Discard the flow-through in the collection tube and return the column into the empty tube.
6. Add 750 µL Buffer PE to the MinElute column and centrifuge at 17,900 RCF for 1 minute; discard flow through and return MinElute column to the collection tube. Centrifuge for 1 minute at 17,900 RCF to remove residual Ethanol.
7. Place MinElute column in a clean 1.5 mL microcentrifuge tube.
8. Add 50 µL EB Buffer (10mM Tris-Cl, pH 8.5) or water directly onto center of the MinElute membrane. Incubate column for 1 minute at RT, and then centrifuge column at 17,900 RCF for 1 minute.
9. Run 2 µL of purified DNA with 5 µL of diluted gel loading dye next to 2 µL pooled unpurified DNA with 5 µL gel loading dye on 0.8% agarose gel for 60 minutes at 100V to verify the purity of the DNA sample with no primer carry-over or contaminating gDNA.
10. This is your purified amplicon pool ready for Illumina sequencing.

### **Suggested Suppliers**

Qiagen: MinElute® PCR Purification Kit (Cat. No. 28004)

Copan: Regular FLOQSwabs® Sterile Collection Swabs 30mm Peel Pouch, Cat #520CS01

~ OR ~

Fisher Scientific: HydraFlock 6" Sterile Elongated Flock Swab w/Polystyrene Handle, 30mm Breakpoint  
Cat No. 22-029-507

Fisher Scientific: Biotech Grade Water BP24854; Potassium Hydroxide P250; TE BP24731

Thermo Fisher: Invitrogen Dyna Mag 96 side magnetic rack, #12331D

Permagen: 0.2 mL PCR Strip Magnetic Separation Rack 8 or 12 Strip, Cat #MSR812

Shoreline Complete™ V4 Protocol 2  
Lyse, Purify, & Amplify Sets C1, C2, C3, C4, C5, C6 (16 tubes) Saliva, Swabs, and Cell Pellets

*This page left intentionally blank*

This document and its contents are proprietary to Shoreline Biome, LLC and are intended solely for the contractual use of its customer in connection with the use of the product(s) described herein and for no other purpose. This document and its contents shall not be used or distributed for any other purpose and/or otherwise communicated, disclosed, or reproduced in any way whatsoever without the prior written consent of Shoreline Biome. Shoreline Biome does not convey any license under its patent, trademark, copyright, or common-law rights nor similar rights of any third parties by this document.

The instructions in this document must be strictly and explicitly followed by qualified and properly trained personnel in order to ensure the proper and safe use of the product(s) described herein. All of the contents of this document must be fully read and understood prior to using such product(s).

**FAILURE TO COMPLETELY READ AND EXPLICITLY FOLLOW ALL OF THE INSTRUCTIONS CONTAINED HEREIN MAY RESULT IN DAMAGE TO THE PRODUCT(S), INJURY TO PERSONS, INCLUDING TO USERS OR OTHERS, AND DAMAGE TO OTHER PROPERTY.**

**SHORELINE BIOME DOES NOT ASSUME ANY LIABILITY ARISING OUT OF THE IMPROPER USE OF THE PRODUCT(S) DESCRIBED HEREIN (INCLUDING PARTS THEREOF OR SOFTWARE) OR ANY USE OF SUCH PRODUCT(S) OUTSIDE THE SCOPE OF THE EXPRESS WRITTEN LICENSES OR PERMISSIONS GRANTED BY SHORELINE BIOME IN CONNECTION WITH CUSTOMER'S ACQUISITION OF SUCH PRODUCT(S).**

All names, logos, and trademarks are the property of their respective owners.